PRO-MLS

[EN] User Guide

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PRO-MLS

Copyright © PROMICRA, s.r.o. 2012


Table of Contents

[EN] PRO-MLS - Universal Heavy Duty Large Stand with Motorized Z Axis	5
1. Preface	5
2. Safety Precautions	6
3. Device Description	7
4. System Assembly	7
4.1. General Assembly	7
4.2. Mounting of Adapters for Optical Devices	8
4.3. Mounting of Microscopic Stages	10
4.4. Mounting of Illuminators	10
4.5. Mounting of Specimen Holders	10
4.6. Cabling	10
5. Limit Switches	11
6. Grounding	11
7. Scheme of Compatibility	12
8. Specification	13
[CZ] PRO-MLS - Univerzální robustní stativ s motorizovanou osou Z určený pro mikroskopické i makroskopické aplikace	15
1. Úvod	15
2. Bezpečnostní pokyny	15
3. Popis zařízení	16
4. Sestavení zařízení	16
4.1. Základní montáž	16
4.2. Montáž adaptérů pro připojení optických zařízení	17
4.3. Montáž mikroskopických stolků	19
4.4. Montáž osvětlovačů	19
4.5. Montáž držáků preparátu	19
4.6. Propojení kabelů	19
5. Koncové spínače	19
6. Uzemnění	20
7. Schéma kompatibility	21
8. Specifikace	22

[EN] PRO-MLS - Universal Heavy Duty Large Stand with Motorized Z Axis

1. Preface

PRO-MLS is an universal motorized stand, which can be used in a wide range of microscopic as well as macroscopic applications. The stand is designed for mounting of a wide range of optical devices as well as microscopic stages and illuminators. The robust construction provides excellent rigidity and high bearing (up to 16kg). The range of Z axis movement is protected by limit switches. The lower limit switch vertical position is adjustable. The stand is designed to be controlled by a CB-ZM-MLS control unit.

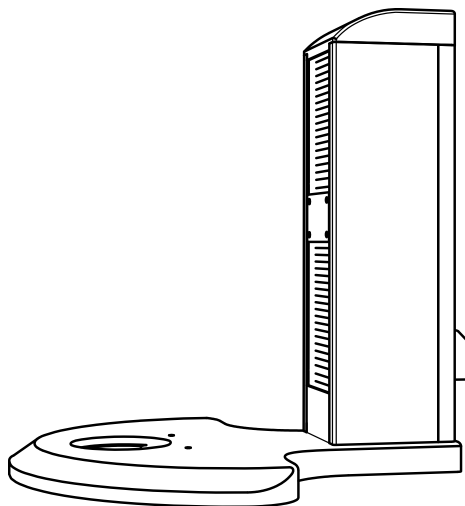2. Safety Precautions

General

1. The moving parts of the PRO-MLS stand are capable of developing a strong force. Never place your hands close to the moving parts of the device. You may take a risk of serious injury.
2. Never open the curtains while the device is in use. Always unplug the cable before entering internal parts of the device.
3. Never stick any objects into the inner spaces of the device while the device is in use.


“CE” mark indicates that this product complies with European EN 61326-1:06 standards.

This symbol (crossed-out wheeled bin WEEE Annex IV) indicates separate collection of waste electrical and electronic equipment in the EU countries. Please do not throw the equipment into domestic refuse. Please use the return and collection systems available in your country for disposal of this product.

RoHS compliance – this product complies with EC directive 2002/95/EC (the RoHS Directive), which restricts the use of the hazardous substances in electrical and electronic equipment.

3. Device Description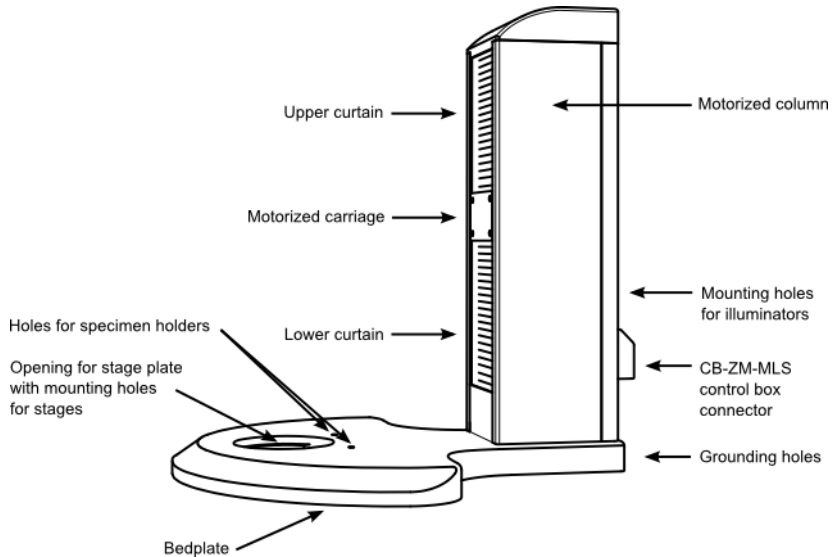


4. System Assembly

4.1. General Assembly

PRO-MLS device is delivered disassembled into two main parts: the motorized column and the bedplate.

The first step of assembly process is the mounting of the motorized column to the bedplate using six M8 hexagon screws (screws and screw driver are included in the package). See figure: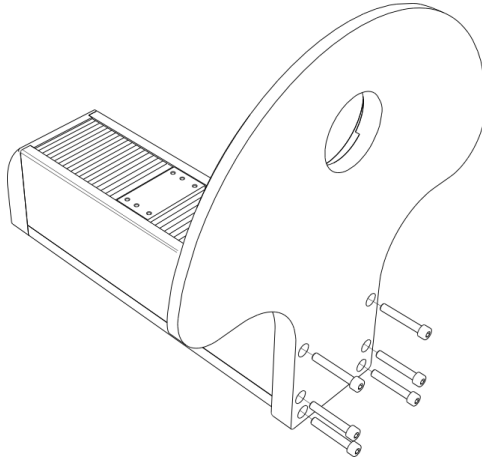

Note

The motorized column does not have to be necessarily mounted to the bedplate. It can be also mounted to a working desk or a table if necessary. This extends rapidly the versatility of the device.


4.2. Mounting of Adapters for Optical Devices

The motorized carriage part of the device is equipped with 6 mounting M5 holes for attaching adapters for various optical devices (microscopes, macro cameras...)

Mounting of PRO-FUAD Adapter


The PRO-FUAD adapter is intended for mounting of OLYMPUS® stereomicroscopes including the focusing unit. For more information see “Scheme of Compatibility“ on page 12.

Mount the PRO-FUAD adapter to the motorized carriage using four hexagon M5 screws and seal the holes by the stickers (screws, stickers and screw driver are included in the adapter package). See figure: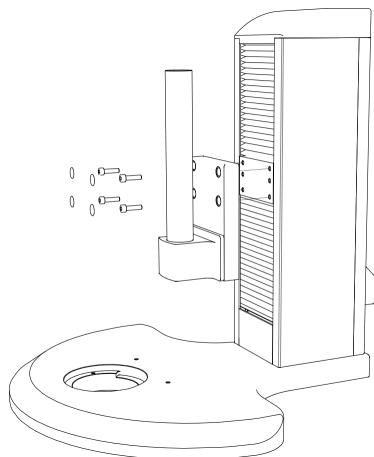

Mounting of PRO-ZBAD Adapter

The PRO-ZBAD adapter is intended for mounting of OLYMPUS® stereomicroscopes without focusing unit. For more information see “Scheme of Compatibility“ on page 12. Mount the PRO-ZBAD adapter to the motorized carriage using four hexagon M5 screws and seal the holes by the stickers (screws and stickers are included in the adapter package). See figure: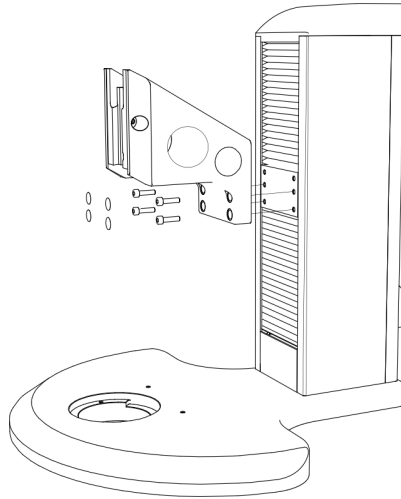

4.3. Mounting of Microscopic Stages

Microscopic stages can be mounted to PRO-MLS stand using two M4 mounting holes located in the stage plate opening.

4.4. Mounting of Illuminators

For mounting of illuminators there are two M4 holes located at the back side of the motorized column.

4.5. Mounting of Specimen Holders

There are two holes located at the top of the bedplate close to the stage plate opening for attachment of specimen holders.

4.6. Cabling

Make sure the CB-ZM-MLS control unit is switched off. Connect the cable included in the package of the CB-ZM-MLS control unit to the connector located at the back side of

the PRO-MLS stand and secure it by two screws. Connect the connector at the second end of the cable to the CB-ZM-MLS control unit and secure it by the screws also.

5. Limit Switches


The PRO-MLS stand is equipped with two limit switches. The limit switches are located inside the motorized column part of the device. The upper limit switch position is fixed. The vertical position of the lower limit switch is adjustable. If there is a microscopic stage mounted to the PRO-MLS stand or specimens of a significant height are to be observed, the lower limit switch position should be adjusted to prevent possible damage of specimens or optics. In order to adjust the vertical position of lower limit switch proceed as follows:

1. Move the motorized carriage to the upper limit using the CB-ZM-MLS control unit.
2. Switch off the CB-ZM-MLS control unit and disconnect the cable from the PRO-MLS stand.
3. Raise up the bottom part of the lower curtain which is fastened by a Velcro fastener. (There is no big force necessary to do this).
4. Locate the limit switch printed circuit board at the left side of the inner space of the motorized column.
5. Loosen up the M6 hexagonal screw which secures the position of the limit switch as well as it works as a mechanical backstop. Do not screw out the screw, just loosen it up.
6. Move the lower limit switch to the desired vertical position.
7. To fix the limit switch position tighten the M6 hexagonal screw.
8. Place the lower curtain to the original position.

6. Grounding

There are four holes located at the rear side of the bedplate which are intended for grounding purposes.

7. Scheme of Compatibility

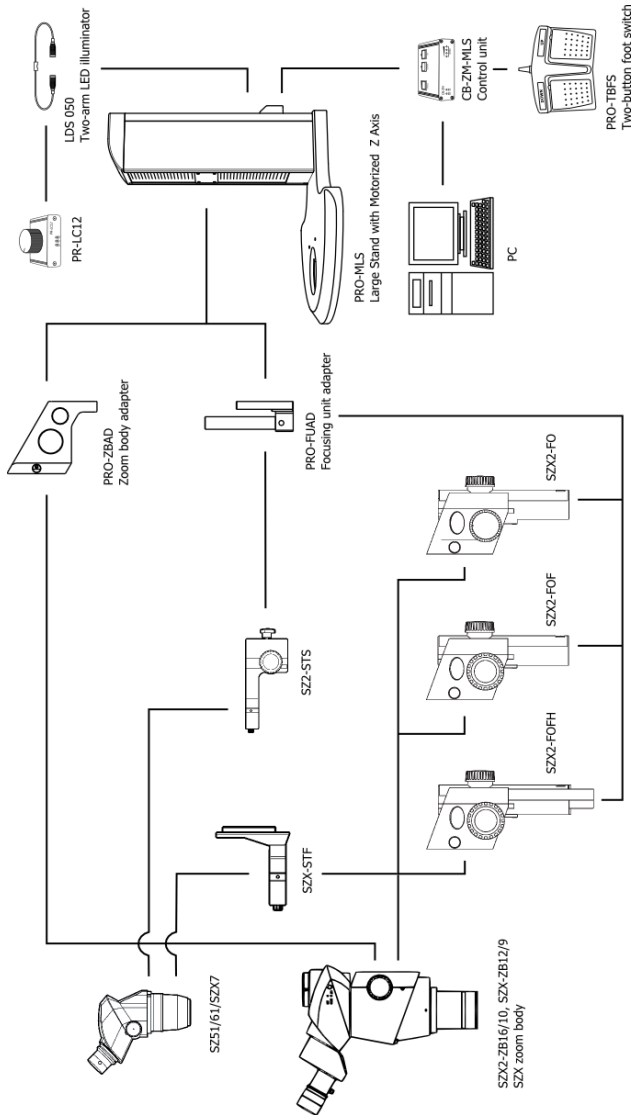8. Specification

Dimensions	45 x45 x47 cm
Control unit	CB-ZM-MLS
Max. bearing	16 kg
Limit switches	lower limit switch adjustable, upper one fixed
Weight	20kg
Manufacturer	PROMICRA, s.r.o., Czech Republic, www.promicra.com

OLYMPUS® is a registered trademark of OLYMPUS Corp.

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OLYMPUS® je registrovaná značka společnosti OLYMPUS Corp.